

Barcode Scanning of Sequel[®] Sequencing Kit 2.0

User Bulletin


PacBio has become aware of an issue with the barcode scanning of the Sequel Sequencing Kit 2.0 (P/N 101-011-000). Scanning of the barcode on the outside of the kit box will allow customers to initially proceed through SMRT[®] Link run set up; however, scanning of the kit barcode will cause a downstream software error that will result in a failed run. Until further notice, customers should scan the barcode on the card located inside the kit instead of the barcode on the box itself.

Procedure

1. Open the Sequel Sequencing Kit 2.0 (P/N 101-011-000).



2. Locate the card containing the barcode.

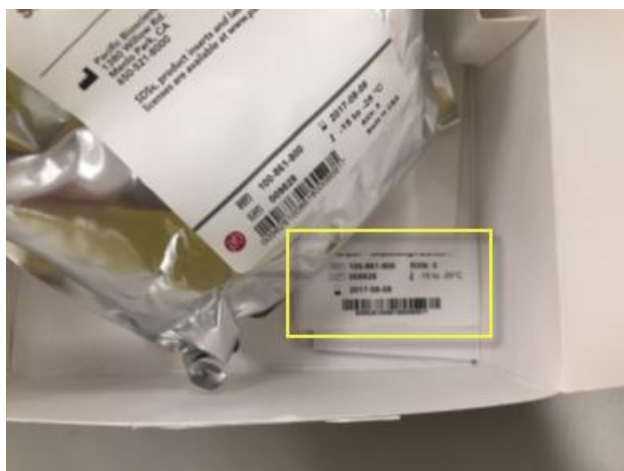

3. Scan the barcode on the card to proceed through SMRT Link.

For Research Use Only. Not for use in diagnostics procedures. © Copyright 2017 by Pacific Biosciences of California, Inc. All rights reserved. Pacific Biosciences, the Pacific Biosciences logo, PacBio, SMRT, SMRTbell, Iso-Seq, and Sequel are trademarks of Pacific Biosciences. BluePippin and SageELF are trademarks of Sage Science. NGS-go and NGSengine are trademarks of GenDx. FEMTO Pulse and Fragment Analyzer are trademarks of Advanced Analytical Technologies. All other trademarks are the sole property of their respective owners.