
Application brief TAXONOMIC AND FUNCTIONAL PROFILING WITH HIFI METAGENOMICS

High-accuracy long-read sequencing to characterize complex microbial communities

HiFi metagenome profiling provides more (and richer) functional information and precise species classifications.


- Each HiFi read typically has an average of eight intact genes per read, providing richer information about genes that co-occur in a single microbe, and may be co-expressed
- 80-90% of HiFi reads have at least one functional annotation, and typically average four, whereas only ~1/3 of short reads can be annotated even once
- · HiFi reads can assign taxonomy with high precision and recall



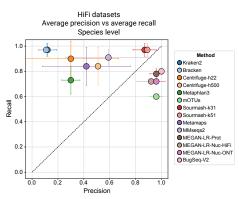

With highly accurate long-read metagenomic data, you can profile species with high precision and recall; tools/methods have been evaluated to provide recommendations to which work best for long-read metagenomic data.¹

Shotgun metagenome functional profiling

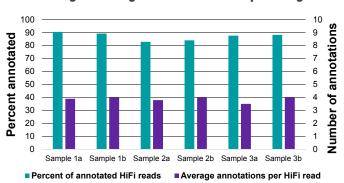

With an average of eight complete genes per HiFi read, PacBio[®] data provides rich functional information; nearly every read contributes to the understanding of the biological functions present in the microbial community.²

Typical workflow from DNA to metagenome taxonomic and functional resolution



Prepare SMRTbell[®] template

Create ~10–15 kb libraries for HiFi metagenome sequencing³



SMRT[®] sequencing and SMRT[®] analysis

Obtain HiFi reads with the Sequel[®] II/e or Revio[™] system



Tertiary data analysis


Profile metagenome taxonomy and functions using our GitHub pipelines⁵ with Sourmash⁶ or Diamond⁷ + MEGAN-LR⁸, or with the BugSeq cloud-based platform⁹

"Our comparisons between long-read sequencing technologies indicate that read quality remains critical for taxonomic profiling performance...The long-read datasets produced significantly better results than the short-read datasets."¹

PacBi

The PacBio partnership with BugSeq empowers public health labs to solve complex microbial questions

Yield top accuracy for taxonomic profiling

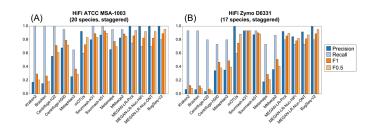

For the best precision/recall trade-off for taxonomic profiling, use BugSeq with HiFi reads.¹ BugSeq offers a curated reference database for analyses requiring the highest quality, and the NCBI nucleotide database for broader taxonomic representation.

Dive deep and analyze antimicrobial resistance

Klebsiella pneumoniae


| .S Copy table III Configure Columns | | I Plot Showin | ng ³¹ / ₃₁ rows and ⁴ / ₄ columns. | | |
|-------------------------------------|----------------------|---------------|--|------------|---|
| Antimicrobial | Class | | Genotypic Predictor of Resistance | Confidence | Genotypic Determina |
| Amikacin | Aminog | lycoside | Present | Very High | aac(6')-lb; aac(6')-lb-cr; aph(3')-VI |
| Streptomycin | Aminog | lycoside | Present | Very High | aadA1 |
| Tobramycin | Aminog | lycoside | Present | Very High | aac(6')-lb; aac(6')-lb-cr |
| Gentamicin | Aminog | lycoside | Not Detected | | |
| Amoxicillin cla | vulanic acid Beta-La | actam | Present | Very High | blaNDM-5 |
| Ampicillin | Beta-La | actam | Present | Very High | blaCTX-M-15; blaNDM-5 blaSHV-11; blaTEM-1A |
| Aztreonam | Beta-La | actam | Present | Very High | blaCTX-M-15 |
| Cefepime Be | | actam | Present | | blaCTX-M-15; blaNDM-5 |
| Cefixime | Beta-La | actam | Present | Very High | blaCTX-M-15; blaNDM-5 |

Demo antimicrobial resistance report from BugSeq outputs. Beyond taxonomic profiling, BugSeq performs *de novo* assembly and identifies predictors of antimicrobial resistance (genes and mutations) as well as virulence factors using curated databases.¹⁰



B 🗏 G S E O

Bioinformatics

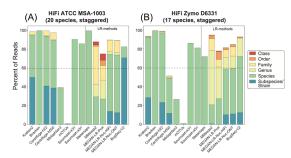

Compatible

In a comparison of analytical tools, BugSeq produced the most accurate abundance estimates and classified the greatest number of reads to the strain rank using PacBio HiFi data.¹

Discover with simplicity

With HiFi data, it's as easy as uploading raw data. BugSeq is the only tool tested by Portik et al. (2022) that is cloud-based, has a graphical user interface, and performs quality control.¹

"Overall, we find that BugSeq, MEGAN-LR-prot, and MEGAN-LR-nuc provide the best tradeoffs for all long-read metagenomics data."¹

- Portik, Brown, and Pierce-Ward, 2022



Learn more about PacBio microbial community applications and BugSeq: pacb.com/metagenomics | bugseq.com/pacbio

KEY REFERENCES

- 1. Portik, D.M., et al. (2022) Evaluation of taxonomic classification and profiling methods for long-read shotgun metagenomic sequencing datasets. *BMC Bioinformatics*, 23: 541. 10.1186/s12859-022-05103-0
- Gehrig, J., et al. (2022) Finding the right fit: evaluation of short-read and long-read sequencing approaches to maximize the utility of clinical microbiome data. Microbial Genomics, 8(3), 10.1099/mgen.0.000794.
- Procedure & checklist Preparing whole genome and metagenome libraries using SMRTbell prep kit 3.0. PacBio documentation.
- Application brochure A new paradigm: HiFi reads for highly accurate long-read sequencing. PacBio documentation.
- https://github.com/PacificBiosciences/pb-metagenomics-tools
- 6. Pierce, N. T., et al. (2019) Large-scale sequence comparisons with sourmash. F1000Research, 8, 1006. 10.12688/f1000research.19675.1
- 7. Buchfink, B., et al. (2015) Fast and sensitive protein alignment using DIAMOND. *Nature methods*, 12(1), 59–60. 10.1038/nmeth.3176
- Huson, D. H., et al. (2018) MEGAN-LR: new algorithms allow accurate binning and easy interactive exploration of metagenomic long reads and contigs. Biology direct, 13(1), 6. 10.1186/s13062-018-0208-7
- https://bugseq.com/ & https://bugseq.com/pacbio
- 10. Chandrakumar et al. (2022) BugSplit enables genome-resolved metagenomics through highly accurate taxonomic binning of metagenomic assemblies. *Communications Biology*, 151(5). 10.1038/s42003-022-03114-4.

Research use only. Not for use in diagnostic procedures. © 2023 Pacific Biosciences of California, Inc. ("PacBio"). All rights reserved. Information in this document is subject to change without notice. PacBio assumes no responsibility for any errors or omissions in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of PacBio products and/or third-party products. Refer to the applicable PacBio terms and conditions of sale and to the applicable license terms at pack.com/license. PacIfic Biosciences, the PacBio logo, PacBio, Circulomics, Omniome, SMRT, SMRTbell, Iso-Seq, Sequel, Nanobind, SBB, Revio, and Onso are trademarks of PacBio.

