



**SMRT Link Chemistry Bundle** updates allow updating of SMRT Link features **without** having to reinstall the SMRT Link software. The update also updates **Sequel® Instrument Control Software (ICS)**.

### Updating the SMRT Link Chemistry Bundle

- **Note:** Only SMRT Link users with the **Admin** role can perform this update.
1. In SMRT Link, choose **About** from the Main menu. (A red circle indicates that a Chemistry Bundle Update is available.)
  2. Click the **Update** button.
  3. Clear the browser cache: Choose **More Tools > Clear browsing data**, choose **the beginning of time** from the droplist, then check **Cached images and files**. Click **Clear browsing data**.
  4. Restart the browser. The SMRT Link Chemistry Bundle for this installation of SMRT Link is updated.
  5. On the instrument, choose **Admin** from the Main menu. (A red circle indicates that a Chemistry Bundle Update is available.)
  6. Click the **Updates** tab, then click **Install**. The instrument software then restarts, which will take around 10 minutes.

### New Features **Sequel® Instrument Control Software (ICS)**

- Enables the use of Sequel Sequencing Kit v2.1 (4 and 8 reactions), which include a surface wetting agent designed to increase SMRT Cell loading performance.

### SMRT Link

- In **Sample Setup**, the on-plate sample loading recommendations were revised.
- **Run Design** now allows selection of the new Sequel Sequencing Kits v2.1.

For Research Use Only. Not for use in diagnostic procedures. © Copyright 2017, Pacific Biosciences of California, Inc. All rights reserved. Information in this document is subject to change without notice. Pacific Biosciences assumes no responsibility for any errors or omissions in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of Pacific Biosciences products and/or third party products. Please refer to the applicable Pacific Biosciences Terms and Conditions of Sale and the applicable license terms at <http://www.pacb.com/legal-and-trademarks/product-license-and-use-restrictions/>.

Pacific Biosciences, the Pacific Biosciences logo, PacBio, SMRT, SMRTbell, Iso-Seq and Sequel are trademarks of Pacific Biosciences. BluePippin and SageELF are trademarks of Sage Science, Inc. NGS-go and NGSengine are trademarks of GenDx. FEMTO Pulse and Fragment Analyzer are trademarks of Advanced Analytical Technologies. All other trademarks are the sole property of their respective owners.