

High-throughput HMW DNA animal blood extraction and sequencing on the PacBio Revio system

Deborah Moine,¹ Adam Bates³, Michelle Kim,² Jacob Brandenburg¹, Nadia Sellami,¹ Jackson Mingle,² Jeffrey Burke,² Julian Rocha,² Aurelie Souppe,¹ Heather Ferrao,¹ Gregory Concepcion,¹ Caroline Howard,³ and Kelvin J Liu.²

1. PacBio, 1305 O'Brien Drive, Menlo Park, CA 94025, 2. PacBio, 701 E. Pratt Street, Baltimore, MD 21202, 3. Sanger Institute, Wellcome Genome Campus, Hinxton, Cambridge, UK

Introduction

PacBie

The Darwin *Tree of Life* project is a large biodiversity initiative aiming to generate high-quality genomes for 70,000 species of eukaryotes across Britain and Ireland. For large-scale projects such as these, high-throughput (HT) solutions are critical; PacBio Nanobind HT DNA extraction kits combined with the Revio system address these needs by significantly increasing throughput and lowering the cost of long-read sequencing.



Samples analyzed

DNA was extracted from the **whole blood** of multiple animals including mammalian and non-mammalian species. For the mammalian animals (bovine, rabbit, porcine), DNA was extracted from 200 μ L blood, whereas for the non-mammalian animals (fish, bird, and reptile), DNA was extracted from 2.5 to 20 μ L.

This difference in input volume is because the red blood cells (RBCs) of mammals do not contain nuclei. In contrast, the RBCs of most other vertebrates are nucleated, thus requiring 10–100× lower blood input volumes in DNA extractions for these animals.

Sequencing on PacBio Revio system

Samples were sequenced on the Revio system at 225 pM to achieve approx. 30× HiFi coverage.

Sample	HiFi yield	Mean HiFi RL	Median QV	Coverage
Bovine 1	82 Gb	14,504 bp	Q35	28×
Bovine 2	106 Gb	15,988 bp	Q34	37×
Minipig	97 Gb	15,750 bp	Q35	36×
Rabbit	93 Gb	15,041 bp	Q35	27×
Chicken 1	98 Gb	16,384 bp	Q34	82×
Chicken 2	91 Gb	16,182 bp	Q34	76×
Tuna	101 Gb	16,759 bp	Q34	128×
Turkey	87 Gb	16,081 bp	Q35	79 ×
Viper	92 Gb	15,414 bp	Q34	65×
Kestrel	83 Gb	16,672 bp	Q34	65×



Workflow

We present fully automated methods for HT high molecular weight (HMW) DNA extraction, shearing, library preparation, and PacBio HiFi sequencing of blood from animals with non-nucleated and nucleated red blood cells (nRBCs).



We can prepare 96 samples from DNA extraction to libraries that are ready for loading in about twelve hours.





Extraction results

Blood samples were extracted in a 96-well plate on Thermo Fisher KingFisher Apex or Hamilton Nimbus Presto systems. Total run time was 2 h 30 min.

Sample name	Taxon	Blood volume	DNA yield	DNA mode size
Bovine 1	Bos taurus	200 µL	15.0 μg	119 kb
Bovine 2	Bos taurus	200 µL	12.2 μg	50 kb
Minipig	Sus scrofa	200 µL	6.8 µg	111 kb
Rabbit	Oryctolagus cuniculus	200 µL	3.9 µg	188 kb
Chicken 1	Gallus domesticus	2.5 µL	11.5 µg	137 kb
Chicken 2	Gallus domesticus	2.5 μL	10.7 μg	87 kb
Tuna	Thunnus orientalis	5 µL	12.5 µg	125 kb
Turkey	Cathartes aura	10 µL	21.0 µg	165 kb
Viper	Squamata viperidae	10 µL	12.0 µg	126 kb
Kestrel	Falconiforms falconidae	20 µL	13.0 µg	102 kb



Rabbit

Mean HiFi

15 Kb







-) HT HMW DNA extraction is performed using a Nanobind CBB HT kit and the ThermoFisher KingFisher APEX or the Hamilton NIMBUS Presto systems. The kit uses Nanobind disks which feature micro- and nano-structured silica wrinkles to shield bound DNA from damage during extraction. Extraction is done on a 96-well plate and can yield from ~4 to 20 µg of HMW DNA in 2.5 hours.
- HT size selection is automated on a Hamilton NGS STAR (protocol under development) to remove DNA <10 kb using the Short Read Eliminator (SRE) kit on unsheared DNA.
- 3) HT shearing + library preparation is also automated by using an automated liquid handler to pipette shear HMW DNA to 15–20 kb (protocol under development). HMW DNA can also be sheared using the Diagenode Megaruptor 3 system. HT library preparation is performed using the SMRTbell prep kit 3.0 (SPK 3.0) on the Hamilton NGS STAR. Manual library preparation can also be performed using the SPK 3.0 kit.

DNA mode size as measured by Femto Pulse is >80 kb for majority of samples — example electropherograms are shown below.





Figure 3. Revio system and HiFi read length distribution plot for 3 samples

De novo assembly

350,000 -

, 300,000

250,000 -

้อ ม_ี 200,000 -

E 150,000 -

For bovine 1 sample, *de novo* assembly was run with 28× coverage obtained with one Revio SMRT Cell using *hifiasm* with default parameter. This bovine assembly shows high contiguity with contig N50 >23 Mb and high quality with accuracy >QV 56. Each haplotype is phased.

	Hap 1	Hap 2
Consensus accuracy	QV 56.4	QV 56.0
N50 contig length	25.4 Mb	23.0 Mb
Max contig length	111.7 Mb	105.5 Mb
Total size	3.04 Gb	2.94 Gb
Number of contigs	727	710

Conclusion

We demonstrated a high-throughput, automated workflow for processing animal blood samples from extraction using Nanobind HT CBB kit through HiFi sequencing on the Revio system.



Figure 1. Nanobind HT CBB kit and Kingfisher Apex automation system (Thermo Fisher)

Figure 2. HMW DNA size distribution on Femto Pulse system (Agilent Technologies)

Library result

were between 15 and 18 kb.

Libraries were prepared following SPK 3.0 workflow automated on Hamilton NGS star or manually. Final libraries DNA mode size measured by Femto Pulse All sequenced samples generated ~30× coverage of HiFi data sufficient for *de novo* assembly.

Acknowledgements

The authors would like to thank Amy Denton for extraction, and Sanger SciOpps team for library preparation and sequencing.

Research use only. Not for use in diagnostic procedures. © 2024 Pacific Biosciences of California, Inc. ("PacBio"). All rights reserved. Information in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of PacBio products and/or third-party products. Refer to the applicable PacBio terms and conditions of sale and to the applicable license terms at pacb.com/license. Pacific Biosciences, the PacBio Solutions of sale and to the applicable PacBio. Circulomics, Omniome, SMRT, SMRTbell, Iso-Seq, Sequel, Nanobind, SBB, Revio, Onso, Apton, and Kinnex are trademarks of PacBio.