

## Preparing Whole Genome Sequencing Libraries using SMRTbell® Prep Kit 3.0 Quick Reference Card

### Additional Information

- [Procedure & Checklist - Preparing Whole Genome Sequencing Libraries using SMRTbell Prep Kit 3.0](#)
- PacBio [Tech Support](#)

### Notes

- This Quick Reference Card requires gDNA inputs ranging from 0.3 µg – 5 µg
- SMRTbell libraries should be stored at 4°C for use within 7 days or -20°C for use after 7 days
- Refer to the Procedure and Checklist for additional protocol details

### Required Materials

SMRTbell Prep Kit 3.0

SMRTbell Cleanup Beads

Low TE Buffer

TempAssure PCR 8-tube strips and compatible magnet

8-channel pipettes for processing multiple samples (200 µL and 20 µL)

Qubit and qubit reagents

8-tube PCR strip tube compatible thermal cycler

1. **Shear 1 µg** of gDNA on the Megaruptor 3 System at a shear speed of **31**.
  - 1.1. Sample should be in EDTA buffer with a volume of **100 – 140 µL** and a target concentration of **20 ng/µL** (acceptable range of 3 ng/µL – 36 ng/µL).
2. Perform bead purification using **SMRTbell Cleanup Beads**.
  - 2.1. **1.0x** volume SMRTbell Cleanup Beads.
  - 2.2. 10-minute incubation at room temperature.
  - 2.3. Wash **2x** with fresh **80% EtOH**.
  - 2.4. Resuspend in **47 µL** of **Low TE buffer** and elute by incubating for **5 minutes** at **room temperature**.
3. Using **1 µL** of sample, dilute and quant the DNA on the Qubit System and measure size on the Femto Pulse System.
4. Perform **Repair and A-tailing** by adding the following to 46 µL of sheared DNA:
  - 4.1. **8 µL** Repair buffer
  - 4.2. **4 µL** End Repair Mix
  - 4.3. **2 µL** DNA Repair Mix
5. Mix well by pipetting and incubate for **30 minutes** at **37°C**, **5 minutes** at **65°C** and **hold** at **4°C**.
6. Perform SMRTbell Adapter Ligation by adding the following reagents in this order to 60 µL of sample from step 5:
  - 6.1. **4 µL** SMRTbell Adapter
  - 6.2. **30 µL** Ligation Mix
  - 6.3. **1 µL** Ligation Enhancer
7. Mix well by pipetting and incubate for **30 minutes** at **20°C** and **hold** at **4°C**.
8. Perform bead purification using **SMRTbell Cleanup Beads**.
  - 8.1. **1x (95 µL)** volume SMRTbell Cleanup Beads.
  - 8.2. 10-minute incubation at room temperature.
  - 8.3. Wash **2x** with fresh **80% EtOH**.
  - 8.4. Resuspend in **40 µL** of EB and elute by incubating for **5 minutes** at **room temperature**.
9. Perform nuclease digestion of incomplete SMRTbell molecules by adding the following to 40 µL of library from step 8.4:
  - 9.1. **5 µL** Nuclease Buffer
  - 9.2. **5 µL** Nuclease Mix

10. **Mix** well by pipetting and incubate for **15 minutes** at **37°C** and **hold** at **4°C**.
11. Immediately proceed to AMPure® PB Beads Size-Selection.
  - 11.1. Make a **35%** dilution of AMPure PB beads by adding **1.75 mL** of resuspended beads to **3.25 mL** of Elution Buffer (can be stored at 4°C for 30 days).
  - 11.2. **3.1x (155 µL)** diluted beads from step 11.1.
  - 11.3. 20-minute incubation at room temperature.
  - 11.4. Wash **2x** with fresh **80% EtOH**.
  - 11.5. Resuspend in **15 µL** of EB and elute by incubating for **5 minutes** at **room temperature**.
12. Using **1 µL** of sample, dilute and quant the DNA on the Qubit System and measure size on the Femto Pulse System.
13. Samples can be stored at **4°C** for short term storage or **-20°C** for long term storage. When
14. ready, proceed to the Sample Setup module in SMRT® Link to prepare SMRTbell libraries for sequencing.

Revision History (Description)	Version	Date
Initial release.	01	Nov 2021

For Research Use Only. Not for use in diagnostic procedures. © Copyright 2021, Pacific Biosciences of California, Inc. All rights reserved. Information in this document is subject to change without notice. Pacific Biosciences assumes no responsibility for any errors or omissions in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of Pacific Biosciences products and/or third party products. Please refer to the applicable Pacific Biosciences Terms and Conditions of Sale and to the applicable license terms at <https://www.pacb.com/legal-and-trademarks/terms-and-conditions-of-sale/>. Pacific Biosciences, the Pacific Biosciences logo, PacBio, SMRT, SMRTbell, Iso-Seq and Sequel are trademarks of Pacific Biosciences. FEMTO Pulse and Fragment Analyzer are trademarks of Agilent Technologies. All other trademarks are the sole property of their respective owners.