

Note: This Instrument Control Software release applies **only** to Revio systems.

Overview

The Revio system Instrument Control Software (Revio ICS) v13.0.0.212033 release is **identical** to Revio ICS v13.0 **except** for the fixed issues noted below. The two versions provide the same run design options and formats, primary and post-primary analysis algorithms, and output file content and formats. We recommend that **all** Revio systems running Revio ICS v13.0 be upgraded to Revio ICS v13.0.0.212033.

Fixed issues

- Fixed an issue that caused the Revio system to operate at lower than the optimal temperature in some circumstances. The lower temperature can slow polymerase activity and so decrease sequencing performance.
- Fixed an issue that causes the Revio system to display “Instrument stage temperature is set out of range” as a false alarm.

Known issues

- At times the touchscreen may freeze and be unresponsive during initialization. The workaround is to hard-reset the instrument to restart initialization by **either** pressing the power button, **or** by unplugging the instrument. Contact PacBio Technical Support if the issue persists.

Note: Contact your FSE (Field Service Engineer) or PacBio Technical Support (support@pacb.com) to perform the Instrument Control Software installation.

Research use only. Not for use in diagnostic procedures. © 2024 Pacific Biosciences of California, Inc. (“PacBio”). All rights reserved. Information in this document is subject to change without notice. PacBio assumes no responsibility for any errors or omissions in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of PacBio products and/or third-party products. Refer to the applicable PacBio terms and conditions of sale and to the applicable license terms at pacb.com/license.

Pacific Biosciences, the PacBio logo, PacBio, Circulomics, Omniome, SMRT, SMRTbell, Iso-Seq, Sequel, Nanobind, SBB, Revio, Onso, Apton, and Kinnex are trademarks of PacBio.

© 2024 PacBio. All rights reserved. For research use only. Not for use in diagnostic procedures.

PN 103-362-100 Version 01 (January 2024)