

This system version supports Sequel II and Sequel Ile systems **only**.

New Features

Support for new sequencing workflows:

- HiFi sequencing now includes 5mC calling in CpG motifs.
- Automatic detection and resolution of heteroduplex inserts in Amplicon applications.
- On-instrument barcode demultiplexing (Sequel Ile system **only**.)
- Adeno-associated virus sequencing mode.

Support for new sequencing consumables:

- SMRTbell[®] prep kit 3.0
- Sequel II binding kit 3.1 and Sequel II binding kit 3.2
- Sequel II sequencing kit 2.0 (1 rxn)
- SMRT[®] Cell 8M single use tray

Reduced data output size: By default, the produced .bam file(s) now contain **only** HiFi reads, resulting in a substantially smaller file size.

Fixed Issues

- Fixed an issue where an alignment camera malfunction could cause a run to abort.
- Fixed an issue with the IUI where certain buttons could be incorrectly disabled.

Research use only. Not for use in diagnostic procedures. © 2022, PacBio. All rights reserved. Information in this document is subject to change without notice. PacBio assumes no responsibility for any errors or omissions in this document. Certain notices, terms, conditions and/or use restrictions may pertain to your use of PacBio products and/or third party products. Refer to the applicable PacBio Terms and Conditions of Sale and the applicable license terms at <http://www.pacb.com/legal-and-trademarks/product-license-and-use-restrictions/>.

Pacific Biosciences, the PacBio logo, PacBio, Circulomics, Omnione, SMRT, SMRTbell, Iso-Seq, Sequel, Nanobind, and SBB are trademarks of Pacific Biosciences of California, Inc (PacBio). All other trademarks are the sole property of their respective owners.